SALE

2017
6/1~7/31

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Reporter Assays · Cell Proliferation/Cytotoxicity Assays · White Microplates

Luciferase Reporter Assays

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
Assay Procedure

britelite plus, neolite, steadylite plus

- 1 Add Reconstitution Buffer to the vial Lyophilized Substrate.

- 2 Add reagent (100 µl/well) to each well.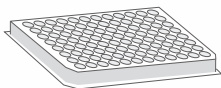


96-well microplate with luc-transformed cells (100 µl/well)

- 3 Mix, wait for 1-15 min. Detect signal.

sensilite

- 1 Add Passive Lysis Solution (20 µl/well) to each well. Shake the plate for 15 min.

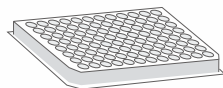
96-well microplate with luc-transformed cells

- 2 Add Reconstitution Buffer to the vial Lyophilized Substrate.


- 3 Inject reagent (100 µl/well) to one well, then detect signal for 5-10 sec. after a 2 sec. count delay. (Repeat this step for next well)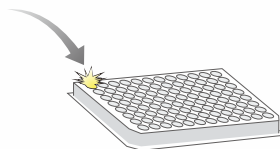

Time course of the luminescent signal

britelite plus shows an intense signal with rapid decrease in signal and neolite and steadylite plus both show strong signal intensity over an extended period of time.

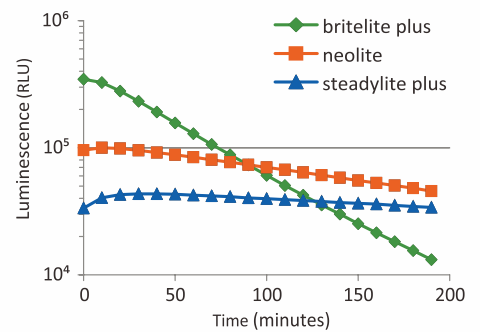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Cell Proliferation and Cytotoxicity Assays

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Luciferase-expression Cell Lines for Drug Screening

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
anti-CD20, anti-HER2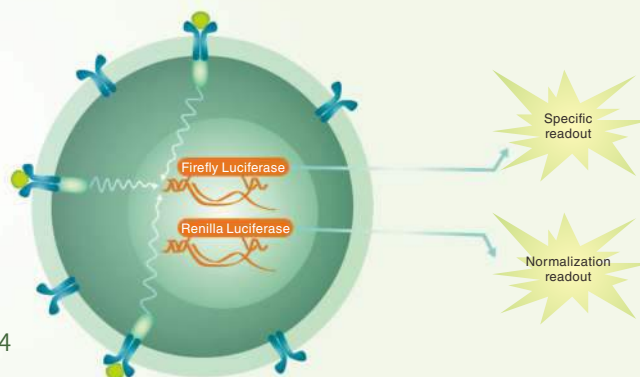

iLite™ cell lines are based on a dual reporter gene technology. The actual drug activity is calculated as the difference between the specific signaling through the target receptor (Firefly Luciferase) and the normalization readout (Renilla Luciferase).